

Algal Toxin Strip Test (Microcystins) 0.3 ppb Finished Drinking Water

1. Collect Sample




Collect 1 to 2 mL of sample.


2. Transfer

Using the pipette provided, add 7 drops of SAMPLE to the conical, flip-top tube.

(The conical, flip-top tube contains dried reagents.)

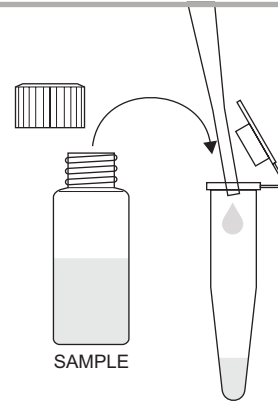

3. Shake and Incubate

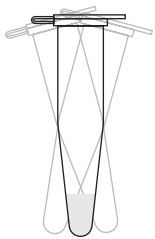

Close the conical, flip-top tube and shake for 30 seconds.

(Dried reagents will dissolve, turning the sample purple.)



Incubate for 10 minutes.



4. Test

Insert test strip into conical, flip-top tube with arrow pointing down.

Incubate for 10 minutes.




5. Dry

Remove test strip. Lay flat and allow to continue developing for 5 minutes.



6. Interpret

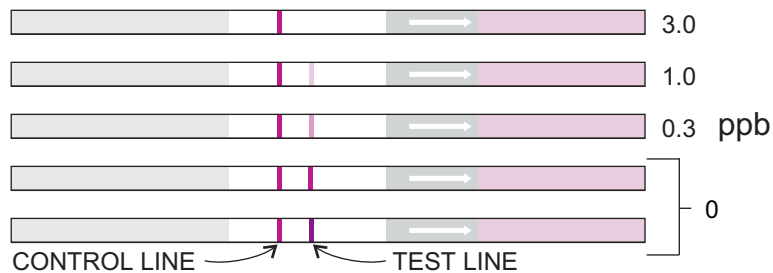

INTERPRET TEST

CONTROL LINE	TEST LINE	INTERPRETATION
NO CONTROL LINE PRESENT	NO TEST LINE PRESENT	INVALID RESULT
CONTROL LINE PRESENT	NO TEST LINE PRESENT	>3 ppb
CONTROL LINE PRESENT	MODERATE INTENSITY TEST LINE PRESENT	BETWEEN 0 AND 3 ppb

For Ordering or Technical Assistance Contact:
 ABRAXIS, INC. 124 Railroad Drive, Warminster, PA 18974
 Phone: 215-357-3911 Fax: 215-357-5232 www.abraxiskits.com

20 Test: PN520046