


Algal Toxin Strip Test (Microcystins) Recreational Water

R110917


1. Collect Sample

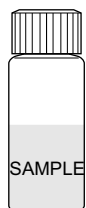

Collect 1 to 2 mL of sample.

2. Transfer/QuikLyse™*

Using the graduated pipette provided, transfer 1 mL of SAMPLE to the lysis vial containing the dried lysis reagent.

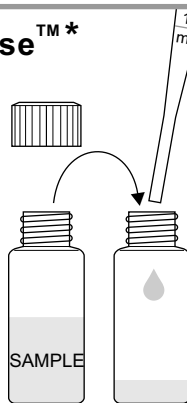

2 min.



8 min.



Cap and shake for 2 minutes.

Let rest for 8 minutes.

*QuikLyse™ reagents may be used in a method of U.S. Patent 9,739,777

3. Add Reagent Paper/QuikLyse™*

Using the forceps provided, add 1 reagent paper to the lysis vial.

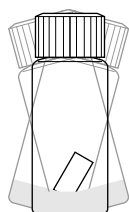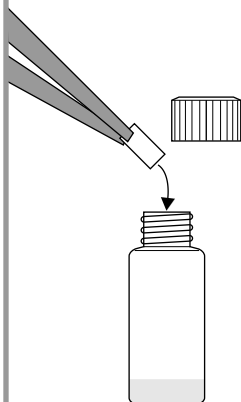

2 min.




8 min.

Cap and shake for 2 minutes.

Let rest for 8 minutes.

4. Transfer


Using the pipette provided, add 7 drops of SAMPLE to the conical, flip-top tube.

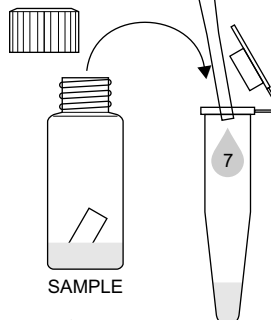

SAMPLE

(The conical, flip-top tube contains dried reagents.)

5. Shake and Incubate

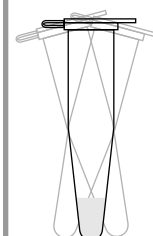

Close the conical, flip-top tube and shake for 30 seconds.



(Dried reagents will dissolve, turning the sample purple.)

6. Test

Insert test strip into conical, flip-top tube with arrow pointing down. (sample pad down).




Incubate for 10 minutes.




7. Dry

Remove test strip. Lay flat and allow to continue developing for 5 minutes.

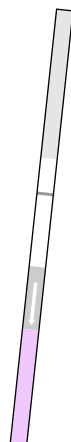

8. Interpret

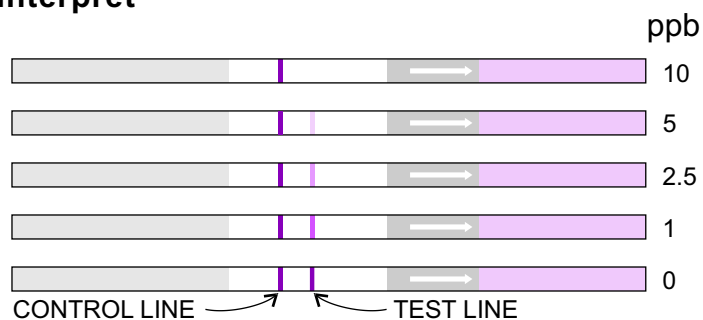

INTERPRET TEST

CONTROL LINE	TEST LINE	INTERPRETATION
NO CONTROL LINE PRESENT	NO TEST LINE PRESENT	INVALID RESULT
CONTROL LINE PRESENT	NO TEST LINE PRESENT	>10 ppb
CONTROL LINE PRESENT	MODERATE INTENSITY TEST LINE PRESENT	BETWEEN 0 AND 10 ppb

For Ordering or Technical Assistance Contact:
ABRAXIS, INC. 124 Railroad Drive, Warminster, PA 18974 Phone: 215-357-3911 Fax: 215-357-5232 www.abraxiskits.com